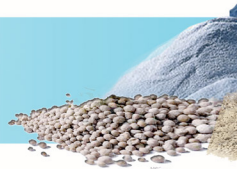
# TUMBLER

*Mixing for Bin*

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*The Tumbler is designed for mixing and homogenization of dry powders, and is primarily used in the Pharmaceutical and Chemical Industries in all those processes of handling of powders in which there is the need to mix homogeneously the powders directly inside Bin without the loading, unloading and cleaning operations of the mixer. The Tumbler produces a well mixing effect thanks to the double 15° inclination of the rotation axis of the machine, and it rotates the BIN both clockwise and counterclockwise.*

*The Tumbler has an electro-mechanical locking system of the bin in order to ensure the horizontal platform position during the loading/unloading of Bin and at the end of each working processing.*

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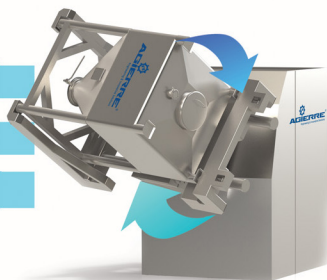
*Automatic closing system with springs and pressure pads system at controlled force for BIN up to 2000 kgs*

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*Mixing and homogenization guaranteed by the double inclination of 15 °*

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*Clockwise and counter-clockwise up to 20rpm*

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*The Tumbler is managed by a PLC system that is connected to the safety sensors and the closing system of the arm. The motor rotates the bin both in a both clockwise and counterclockwise at a speed variable between 4 rpm and 20 rpm.*

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*The wide range of Mixing System shows the great flexibility of our products that can be used in any application within the Pharmaceutical Industry in the handling of Bin or container of any size and shape until 2000 kg a maximum capacity.*

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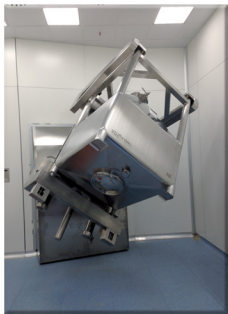
*The Tumbler is designed according to the cGMP standards norms and in accordance with the current safety regulations.*

*It is provided with a set of magnetic sensors and photoelectric barriers that guarantee both the passive and active safety simultaneously.*

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*The Tumbler is designed with a high-tech design and a structure that makes it easy to clean and to look inside.*

*It consists of:*

- **Supporting structure:** it consists of carbon steel profiles anchored to the floor by means of anchor bolts.
- **Lifting platform:** it consists of a stainless steel AISI 304 arm connected to the sliding guides moved by the screw-nut screw system driven by an electric motor.
- **External Carter:** it consists of a stainless steel AISI 304 carter shaped to cover the supporting structure of the Tumbler.
- **Gear motor:** it consists of an electric motor coupled to a reduction gear for the rotation function of the bin.

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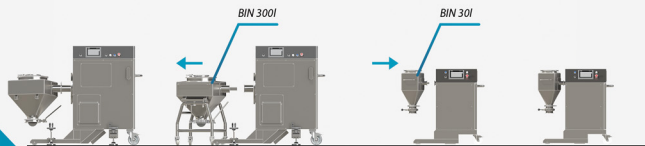
*The fixing and connection system of the Bin can be personalized according to the specific requests of the customer.*

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This type of Bin Blender consists of a Mobile Mixer for BIN designed for mixing and homogenizing dry powders. It is mainly used in the Pharmaceutical and Chemical Industry and in all those handling processes where there is a need to mix the powders homogeneously and directly into the Bin eliminating the loading, unloading and cleaning of the mixer.

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The Bin Blender produces an efficient mixing effect thanks to the 30° inclination of the bin respect to the rotation axis of the machine.

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The Bin Blender is operated by a PLC system, which is connected to the safety sensors and to the locking system of the external structure, which activates the electric motor for the rotation of bin. This motor can rotate the bin clockwise or counterclockwise at a variable speed between 4 rpm and 20 rpm.

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If the Bin Blender is not installed in a working area dedicated exclusively to that machine, it can be equipped with an external structure with an electro-locking system to ensure the safe mixing.

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*The Bin Blender is designed with a highly technological design and a structure that makes it easy to make maintenance activities and clean. It is made of the following components:*

- *Galvanized carbon steel internal structure that represents the machine carrier structure. The arm which support load and the gearmotor are fixed on this structure.*
- *External stainless steel AISI 304 Casing.*
- *Stainless steel AISI 304 security Frame with lexan panels which encompass the bin to be mixed. The entry in the mixing area is allowed by two doors equipped with electro-locking system.*
- *Anti-static wheels with stainless steel AISI 304 support that allow the easy movement of the machine within the department; The rear casters are fitted with a brake.*
- *Stainless steel AISI 304 Arm designed to blend Bin of 5 liters to 300 liters, with a maximum capacity of 150 kgs.*
- *Gearmotor group.*
- *Touch panel from which it's possible to carry out the loading/unloading and mixing of the bin.*

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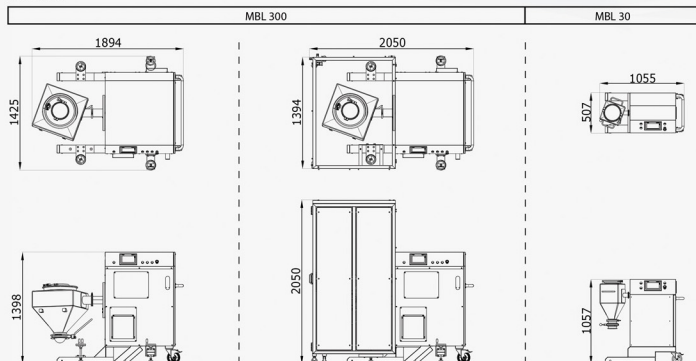
*The Bin Blender is designed in full compliance with current "cGMP" standard norms and in compliance with current safety regulations, UNI EN ISO standards and ATEX regulations (when required).*


## Dimensioni

# Mixing for bin

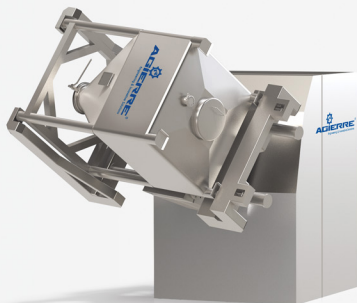
## Tumbler

Per bin fino a 2500l  
For bin up to 2500l

## BL 1500

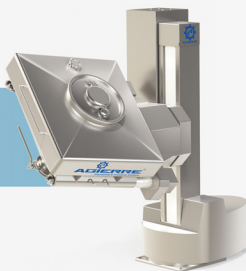Per bin fino a 1500l  
For bin up to 1500l

## MBL 300

Per bin fino a 300l  
For bin up to 300l

## MBL 30

Per bin fino a 30l  
For bin up to 30l